



GenoSensor Corporation

GenoSensor RNA Viral Infection RT-PCR Kit

Catalog # 3401

Version A
February 2024

User's Manual

GenoSensor RNA Viral Infection RT-PCR Kit Manual

Table of Contents

Notes for Instructors	2
Shipping, Storage and Safety	3
GenoSensor RNA Viral Infection RT-PCR Kit Overview	4
Kit Components and Storage Conditions	4
Additional Required Materials	4
Student Guide	6
Full Protocol.....	9
Lab Setting	9
Reagent Preparation	9
RT-PCR Reaction.....	9
Agarose Gel Electrophoresis	10
Results and Discussion	12
Introduction to PCR	14
Troubleshooting	19
Technical Service	20

Literature Citation

When describing a procedure for publication using these products, please refer to them as the *GenoSensor RNA Viral Infection RT-PCR Kit*.

Notes for Instructors

Kit Components and Storage Conditions:

Component	Storage
2X RT-PCR Master Mix	-20°C
Sample A	-20°C
Sample B	-20°C
Sample C	-20°C
Sample D	-20°C
DNA ladder	-20°C

Preparation for RT-PCR (for 6 teams)

1. Set up thermal cycler and the RT-PCR program.
2. Thaw 2X RT-PCR Master Mix on ice.
3. Spin Master Mix for 10 seconds, then vortex for 10 seconds, spin for another 10 seconds.
4. Label 6 microcentrifuge tubes "MM" and aliquot 40 μ L of 2X RT-PCR Master Mix into each tube. KEEP ON ICE.
5. Label 6 tubes (24 total) each "A, B, C, and D" and aliquot 10 μ L of each RNA sample, store on ice.
6. In class, distribute 1 each "MM, A, B, C, D" tubes to all teams.
7. Students will use 10 μ L of 2X RT-PCR Master Mix with 10 μ L sample RNA for a final RT-PCR volume of 20 μ L.

Electrophoresis

- Electrophoresis reagents are not provided in the kit. Please refer to the Additional Required Materials list, on page 4.
- Best results are obtained by adding DNA dye (i.e., Gel Red, Sybr® Safe) to molten agarose.
- For light sensitive DNA dyes, avoid exposing the agarose gel to light. It is best to store and run the gel in a dark room, or cover the gel with a box during gel polymerization and the whole electrophoresis process.
- DNA ladder supplied is enough for 3 lanes with 10 μ L each.
- RNA sample "A, B, C, and D" supplied is enough for up to 20 μ L / lane for 6 teams.

Shipping, Storage and Safety

Shipping and Storage

GenoSensor RNA Viral Infection RT-PCR kits are shipped on blue ice. Components should be stored at temperatures shown in the above table. At proper storage conditions, components are stable for 1 year from the date received. Expiration dates are also noted on product labels.

Safety Warnings and Precautions

This product is intended for research use only. It is not recommended or intended for the diagnosis of disease in humans or animals. Do not use internally or externally in humans or animals. Consider all chemicals as potentially hazardous. Only persons trained in laboratory techniques and familiar with the principles of good laboratory practice should handle these products. Wear suitable protective clothing such as laboratory coats, safety glasses, and gloves. Exercise caution to avoid contact with skin or eyes: if contact should occur, wash immediately with water and follow your laboratory safety protocols. Safety Data Sheets for products are available upon request.

GenoSensor RNA Viral Infection RT-PCR Kit Overview

The GenoSensor RNA Viral Infection RT-PCR Kit introduces common techniques used in RNA research and in medical diagnostics. The kit creates a scenario in which an RNA virus has infected many people and caused a disease. The main symptom of the disease is high fever. Four patients come in a hospital emergency room and they all have high fever. The RNA virus test needs to be done on those patients to diagnose who is infected with the virus. The specific virus gene is going to be tested using this kit and lab results will answer who (A, B, C, and D) has the virus. The test uses reverse transcription polymerase chain reaction (RT-RT-PCR) technology on all four samples. After completing the experiment students will understand the concepts behind RNA detection, RNA amplification, gel electrophoresis, and the genetic concepts driving the experiment.

Kit Components and Storage Conditions

Materials for 6 teams

Component	Amount (27 rxns)	Storage
2X RT-PCR Master Mix	240 μ L	-20°C
Sample A	60 μ L (6 rxns)	-20°C
Sample B	60 μ L (6 rxns)	-20°C
Sample C	60 μ L (6 rxns)	-20°C
Sample D	60 μ L (6 rxns)	-20°C
DNA ladder	30 μ L	-20°C

Additional Required Materials

- Thermal Cycler
- Heat Block or (heat plate, Beaker with de-ionized water; water bath, Tube floater; Thermometer)
- Ice
- Microcentrifuge
- Microcentrifuge tubes (6)
- Vortexer (optional)
- Micropipettes (p10, p200, p1000)
- Pipette tips
- PCR tubes (27)
- Tube Racks
- Ethanol or ethanol wipes
- Electrophoresis equipment
- Microwave oven
- Electrophoresis supplies: agarose, TBE buffer, DNA loading buffer, running buffer, gel dye (e.g., SYBR® safe, Gel Red)

- UV light box or “Gel Doc” equipment and program
- Permanent marker

Student Guide

Objective overview

1. Understand how RNA is responsible for genotypic differences between the test samples of interest.
2. Investigate techniques used in RNA technology: RNA sequence diversity and uniqueness, RNA amplification, RT-RT-PCR, and gel electrophoresis.
3. Investigate and understand the process for gel electrophoresis data analysis.

Many viruses are Ribonucleic acid (RNA) viruses. An RNA virus uses RNA instead of DNA as its genetic materials and can cause many human diseases such as SARS- CoV-2 virus causing the pandemic COVID-19, influenza virus, Ebola virus, HIV (human immunodeficiency virus) and Dengue virus.

RNA is a molecule that is present in the majority of living organisms and viruses. It is made up of nucleotides, which are ribose sugars attached to nitrogenous bases and phosphate groups. The nitrogenous bases include adenine (A), guanine (G), uracil (U), and cytosine (C). RNA mostly exists in the single-stranded form, but there are special RNA viruses that are double-stranded. The RNA molecule can have a variety of lengths and structures.

RNA mutation is a major concern, which can disrupt the normal functioning of RNA and cause potentially life-threatening diseases. The mutation rates of RNA viruses that cause various illnesses in humans are very high. This increase accounts for their fast evolution and ability to produce newer variants with higher infectivity. SARS- CoV-2 that causes pandemic COVID-19 was a good example. Newer strains of SARS- CoV-2 had been identified at a fast pace. The newer variants may also increase resistance to antivirals. Since such mutations are heritable, they provide a bottleneck for the development of drugs or vaccines to combat viral infections. HIV infections are an example of the emergence of many drug-resistant strains, where the viruses replicate and cause severe disease even in the presence of drugs.

RNA detection technologies allow us not only to analyze RNA viruses but also to study many diseases such as cancer diagnosis, diabetic cardiovascular complications, and hematologic malignancies through gene transcript levels and splicing variants. It is important to learn RNA detection technologies.

In this lab you will learn an abridged version of an RNA amplification process. During the exercise you will learn to analyze and compare a number of RNA fragments to determine whether or not the tested RNA fragments are present in the samples. These fragments can be visualized through a process known as "gel electrophoresis".

RNA cannot be directly amplified. RNA needs to be reverse transcribed to its complimentary form of DNA, called cDNA, using reverse transcriptase. Polymerase chain reaction (PCR) can then be applied to amplify this cDNA (Figure 1 and 2). This technology to amplify an RNA is called reverse transcription polymerase chain reaction (RT-PCR).

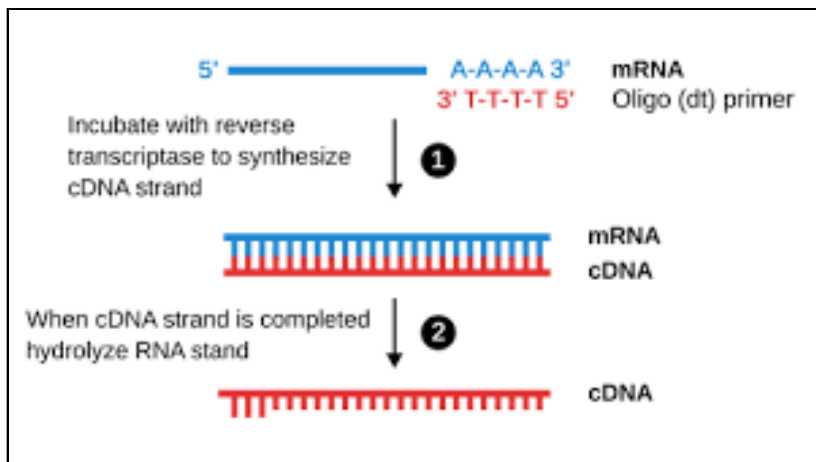


Figure 1, Complementary DNA (cDNA) is reverse transcribed from an RNA by reverse transcriptase.

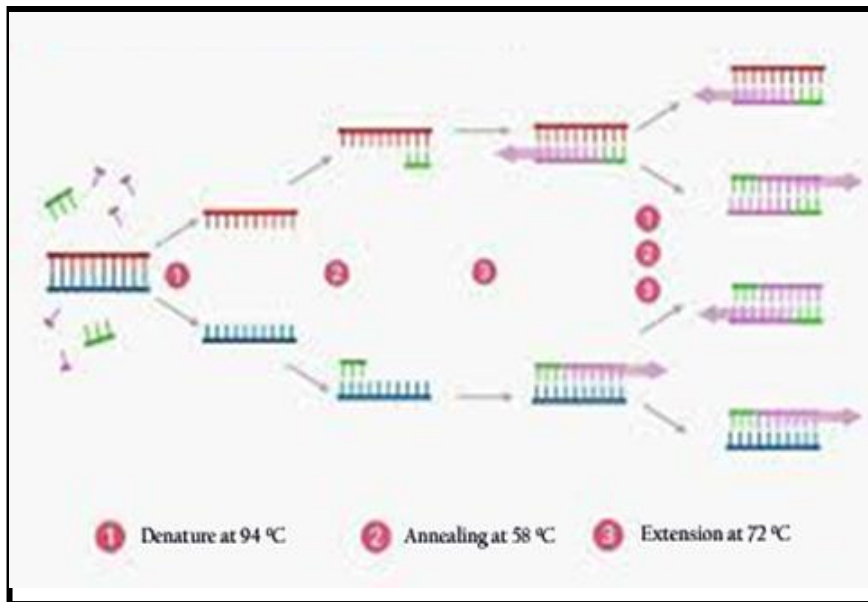


Figure 2, Diagram of PCR process. PCR makes millions to billions of copies of a specific DNA fragment.

RT-RT-PCR is a common technique for life science research, and is also a powerful tool to amplify RNA. Amplification has many purposes and one of them is to see if your sequence of interest is there. If you see amplification of the products, then you can say that the sequence you are interested in is there.

Traditional RT-RT-PCR method uses two steps, reverse transcription and RT-PCR. Reverse transcription step mixes RNA sample with reverse transcriptase and reaction buffer. Incubate the reaction for some time and then take synthesized cDNA for the next RT-PCR step. Nowadays, one-step RT-RT-PCR becomes more popular. RNA sample is mixed with reverse transcriptase, DNA polymerase and optimized buffer together. Incubation of RNA and allowance of reverse transcription take place first. Then thermocycling reaction for the RT-PCR starts after. Reverse transcription occurs at a single and fixed temperature, typically ranging 37 °C – 65 °C.

This kit uses current technology – one-step reaction, and reverse transcription occurs at 46 °C. The amplified products will be examined using gel electrophoresis and UV light. Students will record the results and report who is infected.

Note: Understand the difference between RNA and DNA. DNA is long double helix polymer that uses deoxyribose rings (sugars) and phosphate molecules as support in its backbone. Attached to the backbone are unique sequences of nucleotides which are often referred to as base pairs. There are two different types of nucleotides: purines and pyrimidines. Adenine (A) and Guanine (G) are both purines because they have two rings in their structures. Meanwhile, Thymine (T) and Cytosine (C) are pyrimidines because they have only a single ring in each of their structures. These nucleotides form a bond with their complementary base pair on the other strand of DNA. This is how the double helix structure is formed that resembles a spiral staircase. Each individual will have different sequences of A, T, G, and C in their DNA. There are highly similar and yet unique sequences of DNA that are used to identify humans by looking at the minute differences in their DNA.

Full Protocol

Lab Setting

Materials are enough for 6 groups.

Preparation

1. Set up and program the thermal (see below).
2. Thaw 2x RT-RT-PCR Master Mix on ice.

Reagent Preparation

Refer to "Notes for Instructors – Preparation for RT-PCR" on Page 2.

Pre-Experiment Observations

1. Describe the samples of RNA (physical properties: color, viscosity, etc.). Can you see the RNA?
2. Is there any observable difference between the samples of RNA?
3. Describe the appearance of the 2X RT-RT-PCR Master Mix? Can you see the enzymes?

RT-PCR Reaction

Keep the 2X RT-RT-PCR Master Mix and all samples on ice when not in use.

1. *Wear gloves and handle solutions carefully.* Spin the "MM" tube for 10sec, vortex for 10sec, then spin for another 10sec.
2. **Using a NEW TIP for each sample**, pipette 10 μ L of the 2X RT-PCR Master Mix, (containing reverse transcriptase, Taq DNA polymerase, nucleotides, primers, and RT-PCR reaction buffer) into the sample tubes "A, B, C, and D," already containing 10 μ L of each DNA and the "N" tube if your team received one.

RT-PCR Reaction Mixtures			
RNA Samples		2X RT-PCR Master Mix	Total Reaction Volume
Sample A [A]	10 μ L	10 μ L	20 μ L
Sample B [B]	10 μ L	10 μ L	20 μ L
Sample C [C]	10 μ L	10 μ L	20 μ L
Sample D [D]	10 μ L	10 μ L	20 μ L

3. Pipette up and down carefully to mix well. Tightly cap each tube. Alternatively, mix the components by gently flicking the tubes with your finger. Arrange the tubes in a microcentrifuge and spin for 5 seconds to force all liquid to the bottom of the tubes. (Be sure the tubes are in a BALANCED arrangement in the rotor).

NOTE: If the teacher did not pre-aliquot samples, add the RNA samples THEN the master mix (enzyme) to your microcentrifuge tubes, changing tips each time.

4. Store the samples on ice until they are ready to be loaded into the thermal cycler.

RT-PCR Parameters

Program the thermal cycler as follows:

1. 46°C – 10 minutes
2. 94°C – 30 seconds
3. 94°C denaturing – 20 seconds}
4. 58°C annealing – 20 seconds} **repeat steps 2, 3, & 4 for 35 cycles**
5. 72°C extension – 30 seconds}
6. 72°C – 5 minutes
7. 4°C – finished / hold

STOPPING POINT – For classes with shorter time periods, the RT-PCR samples should be stored at 4°C until the next lab period.

Part 1: Questions

While waiting for the samples, answer the following questions:

1. After combining the 2x RT-PCR master mix with the RNA samples, was there any visible change or any sign of reactivity?
2. Was there any evidence indicating that your samples of RNA were amplified or altered in any way by the addition of the 2X RT-PCR master mix? Explain.

Agarose Gel Electrophoresis

General Procedure, detailed directions as given by instructor

1. Prepare 1% agarose.
2. Set up electrophoresis apparatus and pour in the 1% molten agarose with DNA dye for gelation.
3. For staining, use a DNA dye which is added directly to the molten agarose. For light sensitive dyes, keep the gel in the dark during gelation, either by performing in a dark room or placing a box over the gel.
4. Use at least 10 µL of RT-PCR product to visualize results by electrophoresis on agarose gel. If gel well volume will accommodate more than 10 µL, a higher volume is preferred. Loading dye has already been added to the sample to ensure that the sample will sink to the bottom of the well and properly enter the agarose gel.

Amplified Samples		Loading Volume
Sample A [A]	20 μ L	10 μ L or more
Sample B [B]	20 μ L	10 μ L or more
Sample C [C]	20 μ L	10 μ L or more
Sample D [D]	20 μ L	10 μ L or more
DNA Ladder		10 μ L

- Run at ~100V for ~20 minutes and stop before loading dye has run off gel. Depending on the DNA dye used, caution may need to be taken to reduce exposure of gel to light.
- Visualize under UV light exposure and record the results manually or by photography.

Suggested Gel Setup: Run 3 gels with 10 wells, 2 teams/gel.


Lane 1-4:	Team "1" samples A, B, C, and D up to 20 μ L each.
Lane 5	10 μ L DNA Ladder
Lane 6-9	Team "2" samples A, B, C, and U up to 20 μ L each.

Part 2: Questions

While waiting for your samples to electrophorese, answer these questions with your group members.

- The electrophoresis apparatus creates an electrical field with positive and negative poles at the ends of the gel. DNA molecules are negatively charged. To which electrode pole of the electrophoresis field would you expect DNA to migrate?
- After the DNA samples are loaded into the sample wells, they are "forced" to move through the gel matrix at different speeds. What size fragments (large vs. small) would you expect to move toward the opposite end of the gel and travel a longer distance most quickly? Explain.
- What do you see moving through the gel? (Tricky question).
- Are visualized bands DNA or RNA?

Results and Discussion

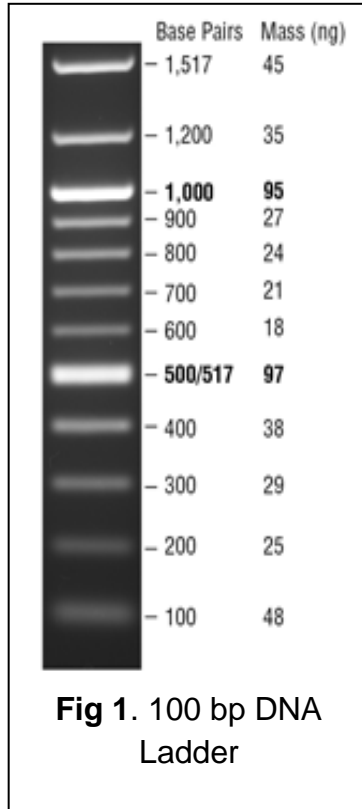

Observe the bands visible in your samples on the gel. Recall which lanes contained the amplified samples and which showed nothing?

Explain why some samples were amplified and some not?

Were all 4 patients infected?

What was happening at the step of 46 °C for 10 minutes?

What was happening at the steps of thermal cycling?

Summarize the process of RT-PCR using the correct terminology.

Describe a new experiment you could perform using RT-PCR and agarose gel electrophoresis.

Final Conclusions

1. Who was infected with the RNA virus?
2. Describe the evidence that confirms the identification of the infectious agent?
3. What are some possible errors that could lead you to have inconclusive results?
4. What further investigations could you perform to better support your evidence?

5. What are the top 3 RNA infectious agents in the US? What diseases do they cause?

6. How else might RT-PCR be used in the medical diagnosis?

Introduction to PCR

In 1983, during his time at Cetus Corporation, Kary Mullis developed a technique that significantly changed the field of genetics and that of all other biological sciences. This revolutionary process was termed “polymerase chain reaction,” or PCR. By 1993, he had earned the Nobel Prize in Chemistry for PCR. His new technique enabled researchers in numerous fields of biology, to easily and rapidly amplify DNA. Before that, amplification of DNA was extremely difficult and time consuming. Now, in the 21st century, it's not just research scientists who use this molecular biology technique. PCR has applications in a wide variety of areas including: gene detection and mapping, whole genome sequencing, analysis of gene expression, forensics, criminal justice, clinical diagnostics, pharmacogenomics, and dozens of others. Nearly every one of these applications was impossible prior to the implementation of PCR. Besides the initial investment in specialized machinery, the cost of performing PCR is relatively low, and the process is simple enough that nearly anyone can do it and get successful results every time!

PCR: The Birth of Recombinant DNA Technology

PCR uses specific nucleotide sequences, named primers, to amplify segments of a genome from a very small amount of starting material, referred to as the template. DNA can be extracted and isolated from almost any cell type i.e., bacterium, blood cells, tissue cells, hair cells, cheek cells, etc. After proper treatment, with PCR, millions of copies of nearly any desired DNA sequence can be produced. The power of PCR is its specificity. PCR uses unique primers to target just the desired sequence of DNA out of the entire genome and amplifies only that segment with little error.

The basic components of PCR:

- Reaction Buffer
- DNA nucleotides (dNTP's) of adenine, guanine, thymine, and cytosine
- DNA polymerase
- Forward and reverse DNA oligonucleotide primers
- Template DNA (starting material)

PCR Makes Use of Two Basic Processes in Molecular Genetics

1. Complementary DNA strand hybridization

For DNA to be amplified, one must have a known sequence that flanks the gene of interest both upstream and downstream. These sequences are used to create ‘oligonucleotide primers,’ meaning a short ~20 base pair nucleotide sequence which is used as a starting point

for DNA replication. The primers are complementary to their target regions so they will anneal (attach) to those regions with great precision. Primers serve the same purpose that runways do for planes trying to lift off the ground. DNA polymerase cannot add nucleotides without a preexisting chain to start from. This process is referred to as primer extension. DNA polymerase recognizes the partially single stranded segment of DNA, attaches itself to the primer just as it normally would during the DNA replication phase of the cell cycle, and proceeds to add complementary nucleotides to fill in the gap.

Complementary strand hybridization occurs when two different oligonucleotide primers anneal to each of their respective complementary base pair sequences on the template. They are designed specifically to anneal at opposite ends of opposite strands of the specific sequence of DNA that is desired to be amplified.

2. DNA strand extension via DNA polymerase

In a PCR, a special type of DNA polymerase is used that is able to function properly and not become denatured during the temperature fluctuation cycles required for thermal cycling. Most mammalian DNA polymerases cannot tolerate the high temperatures and fluctuations from ~60°C-94°C. The breakthrough in PCR came with the isolation of DNA polymerase from a thermophilic bacterium known as *Thermus aquaticus*. This bacterial species lives in high temperature steam vents and its DNA polymerase has evolved to withstand the high temperatures of its environment.

During PCR, DNA is synthesized and its quantity doubles after each cycle making the reaction product grow at an exponential rate. In theory, after 30 cycles there will be 2^{30} -over a billion- copies of DNA. Yielding this much DNA allows it to be visualized after only a few simple procedures. One of the easiest and most popular methods of doing this is agarose gel electrophoresis.

Genes and DNA

The human genome contains 23 pairs of chromosomes that contain a total of thirty to fifty thousand protein coding genes. However, those genes only comprise about 5% of the genome, leaving 95% of it to be classified as non-coding DNA. This non-coding DNA is found not only between, but within genes, splitting them into segments. In eukaryotes, non-coding DNA sequences found within genes are known as **introns**. The sequences that do code for proteins are called **exons**. In eukaryotes, genomic DNA is transcribed into RNA molecules in its unmodified form containing both introns and exons from a particular gene. While the RNA is still in the nucleus (before being transported out of the nucleus), the introns (which **interfere** with the gene product) must be removed from the RNA while the exons (excised from the original transcript) are spliced together to form the complete messenger RNA sequence which will soon be translated into a protein. This process is called RNA splicing. Some genes may contain a few introns, others may contain dozens. Interestingly, it is the non-coding 'junk' DNA that is useful to

us when considering the DNA profile of an individual, rather than the protein coding DNA previously thought to be the most aspect of the genome.

As discussed, functional segments of genes (exons) code for proteins. Proteins are molecules that carry out most cellular functions. Exon sequences are therefore very similar among individuals. That is why when mutations in protein coding genes arise the result is very evident, and is sometime catastrophic. Many metabolic disorders and rare diseases are caused by mutated and nonfunctional proteins). Introns, however, often vary in size and number among individuals. Intron sequences are thought to be the result of the differential accumulation of mutations over time and through evolution are silently passed on to descendants after reproduction. It is differences in the presence and number of intron sequences that allow us to determine the diversity of human genetics. The recognition of these distinctive characteristics in DNA represents the molecular basis from which human identification and population genetics are made possible. Throughout evolution, intron sequences have been the target of random insertions by short repetitive interspersed elements (SINEs), or long repetitive interspersed elements (LINEs). Those elements have become randomly inserted within our introns over millions of years.

PCR Stages

The machinery required to perform PCR is known as a thermal cycler. The thermal cycler enables the steps of PCR to be automated. The reaction involves a repetition of cycles that promote template denaturation, primer annealing, and primer extension by the Taq DNA polymerase. A DNA sample is added to a mixture of the necessary reagents: oligonucleotide primers, thermostable DNA polymerase (Taq), the four nucleotides (A, T, G, C), and reaction buffer. These reagents are pre-mixed as a 2X PCR Master Mix in the GenoSensor Food Poisoning kit II. The tubes are placed into the thermal cycler which contains an aluminum block that holds the samples and can be rapidly heated and cooled by extreme temperature changes in a controlled environment. The rapid heating and cooling of this thermal block is called temperature cycling or thermal cycling.

The first step of the PCR temperature cycling procedure heats the sample to 94°C causing the template strands separate. This process is called **denaturation**.

The thermal cycler then rapidly cools to 58°C allowing the primers to anneal to the separated template strands. This is the **annealing** process. The two original template strands may re-anneal to each other or compete with the primers at the primers' complementary binding sites. However, the primers are added in excess so that the primers may out-compete the original DNA strands for their complementary binding sites.

Lastly, the thermal cycler heats the sample to 72°C (the usual environment temperature for *Thermus aquaticus*) so that Taq DNA polymerase can perform **primer extension** and produce complementary DNA strands of the target sequence. The two resulting new sets of

double-stranded DNA (dsDNA) will be used for the next cycle and proceeding strand synthesis. At this stage, a full temperature cycle (thermal cycle) will have been completed.

Each step takes 30 seconds to 1 minute, and will repeat for 30-40 cycles depending on how the user has programmed the thermal cycler. At the end of the pre-programmed number of cycles, the product is put on hold at 4°C until the user is ready to retrieve the PCR product and analyze its contents.

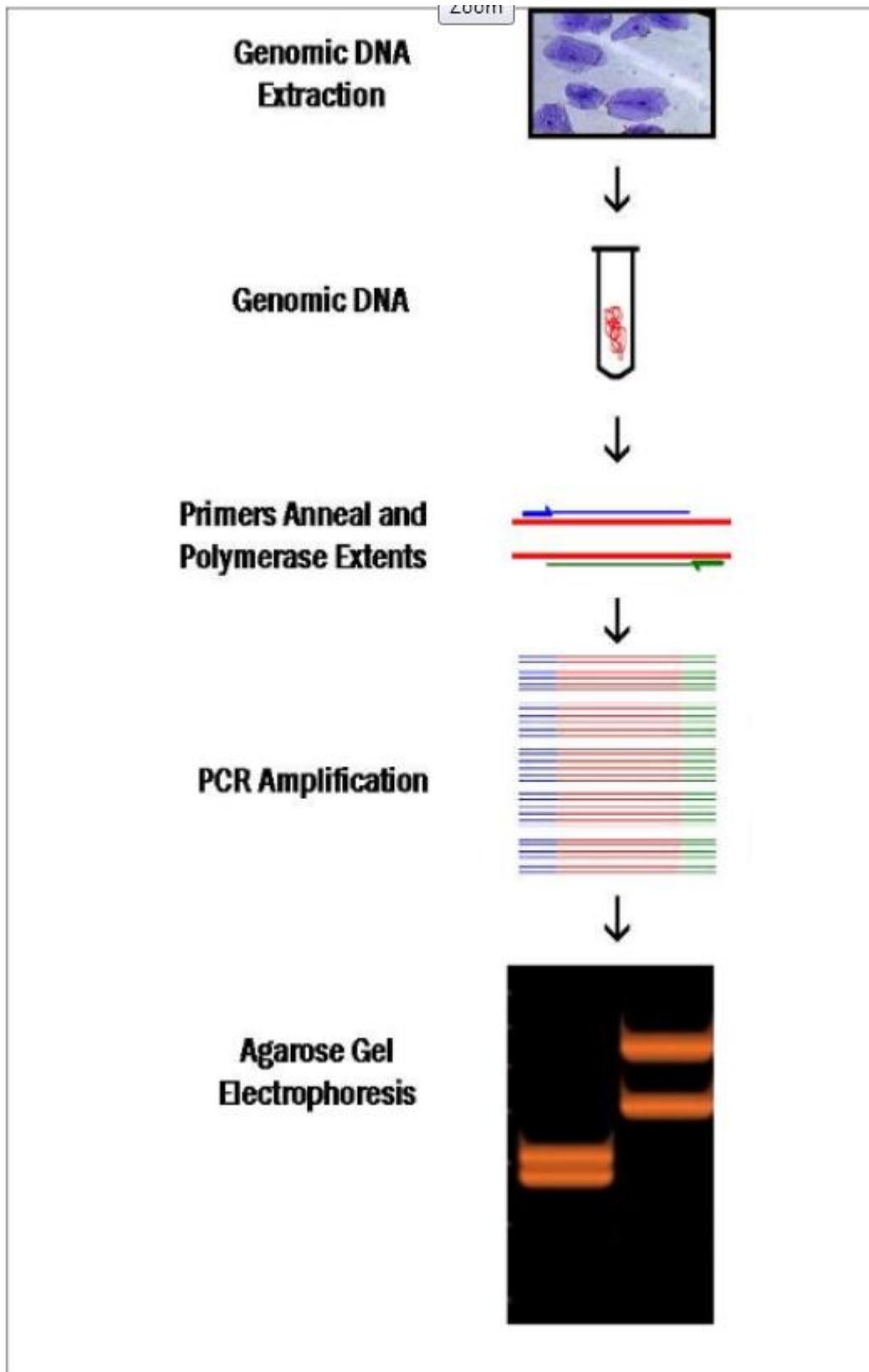


Figure 3. Experiment flowchart from start to finish

Troubleshooting

Symptom	Possible causes	Solutions
No amplification product	Questionable template quality	Analyze starting material
	Inhibitory Substance in reaction	Decrease sample volume
	Insufficient cycle #	Run additional cycles
	Incorrect thermal cycler program	Verify times and temperatures
	Errors in heat block incubation	Calibrate heating block, use sand or water to maximize contact with tube for proper heat transfer
	Contaminated tubes/solutions	Autoclave tubes and use filter tips
	Primer annealing temperature too high	Lower annealing temperature in 2°C increments
Weak bands/faint signal	Low concentration of DNA template	Make sure enough DNA has been added; see recommended amount of DNA to load into RT-PCR reaction.
	DNA Dye degradation during preparation	Light sensitive dyes should be kept in the dark during gel preparation. Prepare in dark room or place a box over the electrophoresis apparatus during gelation and electrophoresis.
	Expired, contaminated or degraded DNA dye	Verify that the DNA dye has not degraded in storage, been contaminated or expired.
	Primer annealing temperature too low	Raise annealing temperature in 2°C increments
	Insufficient mixing of reaction solution	Mix solutions thoroughly before beginning the reaction
	Exogenous DNA contamination	-Wear gloves -Use dedicated area for sample preparation -Use non-aerosol tips

Technical Service

For more information or technical assistance, please email.

GenoSensor Corporation
4665 S. Ash Avenue
Suite G-18
Tempe, Arizona 85282
Email: tech_service@genosensorcorp.com
Web: www.genosensoreducation.com

Limited Warranty

GenoSensor is committed to providing our customers with high-quality goods and services. Our goal is to ensure that every customer is 100% satisfied with our products and our service. If you should have any questions or concerns about a GenoSensor product or service, please contact our Technical Service at tech_service@genosensorcorp.com. GenoSensor warrants that all of its products will perform according to the specifications stated on the certificate of analysis. This warranty limits GenoSensor Corporation's liability only to the cost of the product. No warranty is granted for products beyond their listed expiration date. No warranty is applicable unless all product components are stored in accordance with instructions. GenoSensor reserves the right to select the method(s) used to analyze a product unless GenoSensor agrees to a specified method in writing prior to acceptance of the order. GenoSensor makes every effort to ensure the accuracy of its publications, but realizes that the occasional typographical or other error is inevitable. Therefore GenoSensor makes no warranty of any kind regarding the contents of any publications or documentation. If you discover an error in any of our publications, please report it to our Technical Service. GenoSensor assumes no responsibility or liability for any special, incidental, indirect or consequential loss or damage whatsoever. The above limited warranty is sole and exclusive. No other warranty is made, whether expressed or implied, including any warranty of merchantability or fitness for a particular purpose.